



# Bovine IgG Assay

## Immunoenzymetric Assay for the Measurement of Bovine IgG Catalog # F070

### Intended Use

This kit is intended for use in quantitating bovine IgG. The kit is for Research and Manufacturing use only and is not intended for diagnostic use in humans or animals.

to the concentration of bovine IgG present in the sample. Accurate quantitation is achieved by comparing the signal of unknowns to bovine IgG standards assayed at the same time.

### Summary and Explanation Of the Test

The manufacture of products by various biotechnological processes such as cell or tissue culture can result in residual contamination of the desired product by components used in the culture media. Many formulations of media contain either bovine serum or other more purified fractions of serum proteins such as albumin or transferrin which may contain varying amounts of bovine IgG. When the intended product may be used as a therapeutic agent in humans or animals the product should be highly purified to avoid potential health risks or other problems that might result from trace contaminants. Efforts to reduce trace media contamination to the lowest level practical through optimal process design, validation, and final product testing require a highly sensitive and reliable analytical method. The *Cygnus Technologies* Bovine IgG assay is designed to provide a simple to use, precise, and highly sensitive method to detect contamination to less than 250pg/mL. As such, this kit can be used as a tool to aid in optimal purification process development and in routine quality control of in-process streams as well as final product.

### Reagents & Materials Provided

Component	Product #
<b>Anti-bovine IgG labeled with HRP</b> Sheep antibody (H&L chain specific) conjugated to horseradish peroxidase with preservative, 1x12mL	<b>F071</b>
<b>Anti-bovine IgG coated microtiter strips</b> 12x8 well strips in a bag with desiccant	<b>F072</b>
<b>Bovine IgG Standards</b> Bovine IgG in a protein matrix with preservative. Standards at 0, 0.25, 1, 4, and 20 ng/mL. 1mL/vial	<b>F073</b>
<b>TMB Substrate</b> 3,3',5,5'-Tetramethyl benzidine, 1x12mL	<b>F005</b>
<b>Stop Solution</b> 0.5N Sulfuric Acid, 1x12mL	<b>F006</b>
<b>Wash Concentrate (20X)</b> Tris buffered saline with preservative, 1x50mL	<b>F004</b>

### Storage & Stability

- \* All reagents should be stored at 2°C to 8°C until the expiration date printed.
- \* The substrate reagent should not be used if its' absorbance (stopped) at 450nm is greater than 0.1.
- \* Reconstituted wash solution is stable until the expiration date of the kit.

### Principle of the Procedure

The Bovine IgG assay is a two-site immunoenzymetric assay. Samples containing Bovine IgG are reacted in microtiter strips coated with an affinity purified capture antibody. A second HRP labeled anti-Bovine IgG antibody is reacted, forming a sandwich complex of solid phase antibody-Bovine IgG-HRP labeled antibody. After a wash step removes any unbound reactants, the strips are then reacted with TMB substrate. Followed by the addition of a stop solution changing the color from blue to yellow. The amount of hydrolyzed substrate is read on a microtiter plate reader and will be directly proportional

### Materials & Equipment Required But Not Provided

- Microtiter plate reader spectrophotometer with dual wavelength capability at 450 & 650nm.  
*(If your plate reader does not provide dual wavelength analysis you may read at just the 450nm wavelength.)*
- Pipettors - 50µL and 100µL
- Repeating or multichannel pipettor - 100µL
- Microtiter plate rotator (150 - 200 rpm)
- Sample Diluent, (recommend *Cygnus* Cat. # F031A)
- Distilled water
- 1 liter wash bottle for diluted wash solution

## Precautions

For research or manufacturing use only.

Stop Reagent is 0.5N H<sub>2</sub>SO<sub>4</sub>. Avoid contact with skin and eyes. Flush exposed areas immediately with water.

## Preparation of Reagents

- \* Bring all reagents to room temperature.
- \* Dilute wash concentrate to 1 liter in distilled water, and label with kit lot and expiration date. Store at 4°C.

## Procedural Notes

1. Complete washing of the plates to remove excess unreacted reagents is essential to good assay reproducibility and sensitivity. We advise against the use of automated or other manual operated vacuum aspiration devices for washing plates as these may result in lower specific absorbances, higher non-specific absorbance, and more variable precision. The manual wash procedure described below generally provides lower backgrounds, higher specific absorbance, and better precision. If duplicate CVs are poor or if the absorbance of the “0” standard minus a substrate blank is greater than 0.2, evaluate plate washing procedure for proper performance.

2. This kit is a very sensitive assay for bIgG (< 150pg/mL). Since bovine serum and other bovine protein products are common reagents in many laboratories and are often used at relatively high concentrations it is very important to use extreme care to avoid contamination of any of the components in this kit with external sources of bIgG or other bovine protein products. bIgG contamination will manifest itself as either high assay background, poor precision, or unexpected results.

3. Dilution of samples will be required for samples >20ng/mL. The diluent used should be compatible with accurate recovery. The preferred diluent is our Cat. # F031A available in 100mL, 500mL, or 1 liter bottles. This is the same material used to prepare the kit standards. As the sample is diluted in F031A its matrix begins to approach that of the standards thus reducing any inaccuracies caused by dilutional artifacts. Other prospective diluents should be validated in the assay to demonstrate that they do not give elevated background and are not contaminated with bIgG or other bovine proteins. The diluent should also give acceptable recovery when spiked with known quantities of bIgG.

4. High Dose Hook Effect may be observed in samples with very high concentrations of bIgG. Samples greater than 2µg/mL may give absorbances less than the 20ng/mL standard. If a hook effect is possible samples should also be assayed diluted. If the

dilution corrected concentration of the diluted sample is greater than the undiluted sample this may be indicative of the hook effect.

## Limitations

- \* The antibodies used in this kit may cross react with bovine IgM. Cross reactivity with other species has not been extensively investigated.
- \* Certain sample matrices may interfere in this assay. Although the assay is designed to minimize matrix interference, materials such as detergents in high concentration, extremes of pH (<6.0 and >8.5) or very high protein concentrations may give erroneous results. It is recommended to test the sample matrix for interference by diluting the 20ng/mL standard 1 part to 3 parts of the matrix that does not contain any bIgG. This diluted standard when assayed as an unknown should give a value of 5ng/mL ± 20%. In cases where bIgG levels in the sample will allow for sample dilution, such dilution will often overcome sample matrix interference. Consult *Cygnus Technologies* Technical Service Department for advice on how to quantitate the assay in problematic matrices.

## Assay Protocol Notes

\*The assay is very robust such that assay variables like incubation times, sample size, and other sequential incubation schemes can be altered to manipulate assay performance for more sensitivity, increased upper analytical range, or reduced sample matrix interference. Before modifying the protocol from what is recommended, users are advised to contact our technical services for input on the best way to achieve your desired goals.

\* The protocols specify use of an approved microtiter plate shaker for the immunological step. If you do not have such a device, it is possible to incubate the plate without shaking however it will be necessary to extend the immunological incubation step in the plate by about 60 minutes in order to achieve comparable results to the shaking protocol. **Do not shake during the 30 minute substrate incubation step as this may result in higher backgrounds and worse precision.**

\* Avoid the assay of samples containing Sodium Azide, (NaN<sub>3</sub>) which will destroy the HRP activity of the conjugate and could result in the under-estimation of bIgG levels in that sample.

\* Bring all reagents to room temperature.

\* Set-up plate spectrophotometer to read dual wavelength at 450nm for the test wavelength, and 650nm for the reference wavelength. Blank the instrument using the zero standard wells after assay completion.

\* All standards, controls and samples should be assayed in duplicate. Samples that could contain very high levels of bIgG above the 20ng/mL standard or in the “Hook” region of this assay should also be assayed diluted. Recommended diluent is *Cygnus Technologies*

Cat # F031A. Avoid the use of diluents which contain  $\text{NaN}_3$  or could be contaminated with trace levels of bIgG or other bovine proteins.

\* Thorough washing is essential to proper performance of this assay. Automated plate washing systems or other vacuum aspiration devices are not recommended. The manual method described in the assay protocol is preferred for best precision, sensitivity and accuracy. A more detailed discussion of this procedure can be obtained from our Technical Services Department or at our web site.

\* Maintain a repetitive timing sequence from well to well for all assay steps to insure that all incubation times are the same for each well.

\* Make a work list for each assay to identify the location of each standard control and sample.

\* If the substrate has a distinct blue color prior to the assay it may have been contaminated. If this appears to be the case, read 100 $\mu\text{L}$  of substrate plus 100 $\mu\text{L}$  of Stop Solution against a water blank. If the absorbance is greater than 0.1 it may be necessary to obtain new substrate or the sensitivity of the assay may be compromised.

\* Plates should be read within 30 min. after adding stop since color will fade over time.

### Protocol

**1. Pipette 50 $\mu\text{L}$  of standards, controls and samples into wells indicated on work list.**

**2. Pipette 100 $\mu\text{L}$  of anti-bovine IgG:HRP (#F071).**

**3. Cover and transfer to rotator. Incubate at 180rpm for 2 hours at ambient temperature, 24°C  $\pm$  4°.**

**4. Dump contents of wells into waste or gently aspirate with a pipettor. Blot and vigorously bang out residual liquid over absorbent paper. Fill wells generously with diluted wash solution by flooding well from a squirt bottle or by pipeting in ~350  $\mu\text{L}$ . Dump and bang again. Repeat for a total of 4 washes.**

**5. Pipette 100 $\mu\text{L}$  of TMB substrate (#F005).**

**6. Incubate for 30 minutes. Do Not Shake!**

**7. Pipette 100 $\mu\text{L}$  of Stop Solution (#F006).**

**8. Read absorbance at 450/650nm blanking on the Zero standard.**

## Example Data

Well #	Contents	Abs. at 450nm	Mean Abs.	ng/mL bIgG
1A	Zero Std	0.000		
1B	Zero Std	0.002	0.001	
1C	0.25ng/mL	0.035		
1D	0.25ng/mL	0.031	0.033	
1E	1ng/mL	0.124		
1F	1ng/mL	0.123	0.124	
1G	4ng/mL	0.467		
1H	4ng/mL	0.485	0.476	
2A	20ng/mL	1.785		
2B	20ng/mL	1.705	1.745	
2C	sample 1	0.005		
2D	sample 1	0.010	0.008	<0.1ng
2E	sample 2	0.127		
2F	sample 2	0.123	0.125	1ng

## Calculation of Results

The standards may be used to construct a standard curve with values reported in ng/mL. This data reduction may be performed through computer methods using curve fitting routines such as point to point, spline, or 4 parameter logistic fit. **Do not use linear regression analysis to interpolate values for samples as this may lead to significant inaccuracies!** Data may also be manually reduced by plotting the absorbance values of the standard on the y-axis versus concentration on the x-axis and drawing a smooth point to point line. Absorbances of samples are then interpolated from this standard curve.

## Quality Control

-Precision on duplicate samples should yield average % coefficients of variation of less than 10% for samples greater than 1ng/mL. CVs for samples <1ng/mL may be greater than 10%.

- For optimal performance the absorbance of the substrate when blanked against water should be <0.1.

-It is recommended that each laboratory assay appropriate quality control samples in each run to insure that all reagents and procedures are correct.

## Performance Characteristics

*Cygnus Technologies* has validated this assay by conventional criteria as indicated below. A more detailed copy of this "Validation Summary" report can be obtained by request. This validation is generic in nature and is intended to supplement but not replace certain user and product specific validation that should be performed by each laboratory. At a minimum each

laboratory is urged to perform a spike and recovery study in their sample types. In addition, any of your samples types containing bIgG within or above the analytical range of this assay should be evaluated for dilutional linearity to insure that the assay is accurate and has sufficient antibody excess for your particular bIgG samples. Each laboratory and technician should also demonstrate competency in the assay by performing a precision study similar to that described below. A more detailed discussion of recommended user validation protocols can be obtained by contacting our Technical Services Department or at our web site.

**Sensitivity**

The lower limit of detection (LOD) is defined as that concentration corresponding to a signal two standard deviations above the mean of 10 replicates of the zero standard. LOD is ~100 pg/mL in the recommended protocol. The lower limit of quantitation (LOQ) is defined as the lowest concentration, where concentration coefficients of variation (CVs) are <20%. The LOQ is ~150 pg/mL.

**Precision**

Precision is defined as the percent coefficient of variation (%CV). This is calculated by dividing the standard deviation by the mean value for a number of replicate determinations of two different control samples in the low and high concentration range of the assay. Both intra and inter-assay (n=5 assays) precision were determined on 2 pools with low (~1ng/mL) and high concentrations (~5ng/mL).

Intra-assay			Inter-assay		
# of tests	Mean ng/mL	%CV	# of assays	Mean ng/mL	%CV
20	1.1	7.2	5	1.1	7.8
20	5.2	4.3	5	5.0	6.6

**Specificity/Cross-Reactivity**

In sandwich ELISA cross reactivity can manifest itself either as a false increase in bIgG (positive cross reactivity) or as a false decrease in bIgG (negative cross reactivity) when bIgG present in the sample competes with the cross reactant for the kit antibodies. Mouse IgG and human IgG1, IgG2a, IgG2b, and IgA were tested at 2 mg/mL and found to have no cross reactivity in this assay. All matrices should be evaluated by the user for the possibility of cross reactivity.

**Recovery/ Interference Studies**

Various buffer matrices were evaluated by adding known amounts of bIgG. Because this assay is designed to minimize matrix interference most of these buffers yielded acceptable recovery defined as between 80-120%. In general, extremes in pH (<6.0 and >8.5), high salt concentrations, as well as certain detergents can cause under-recovery. Some product proteins in high concentration may also interfere in the accurate

measurement of bIgG. Each user should validate that their sample matrices yield accurate recovery. Such an experiment can be easily performed by diluting the 20ng/mL standard provided with this kit into the sample matrix in question. For example, add 1 part of the 20ng/mL standard to 3 parts of the matrix containing no or very low bIgG contaminants. This diluted standard when assayed as an unknown should give a value of 5 +/- 20%. Consult *Cygnus Technologies* Technical Service Department for advice on how to quantitate the assay in problematic matrices.

**Hook Capacity**

Increasing concentrations of HCPs > 20 ng/mL were assayed as unknowns. The hook capacity, defined as that concentration that can give an absorbance reading less than the 20ng/mL standard was 2 µg/mL.

**Ordering Information/  
Customer Service**

**To place an order or to obtain additional product information contact *Cygnus Technologies* Customer Support:**

**Tel: 910-454-9442 Fax: 910-454-9443**  
**Email: cygnustec@aol.com**  
**Web: www.cygnustechnologies.com**  
**4701 Southport Supply Road SE, Suite 7**  
**Southport, NC 28461 USA**